



- **NOCOBX**® are Petri dishes with nutrients aimed to highlight microorganism presence on surfaces. It is used to check the result obtained by automated surface disinfection devices.
- **NOCOBX**® contain nutritive substance that promotes the growth of bacteria and therefore highlight their presence.

REFERENCES AND PACKAGING

Reference	Packaging
5000.020	Box of 10 Petri dishes

STORAGE

- Store the product in a cool (temperature below 30°C) and dry place.
- Avoid exposing boxes to sunlight.

INSTRUCTIONS FOR USE

- Open the box carefully (do not touch the test area).
- Apply 10 seconds with a light pressure against the surface to be tested.
- Close the box.
- Leave the box at development temperature, 20-25°C for 48 hours.
- Take the reading.

IMPORTANT :

- Elimination: treatment and incineration by specialized companies.

OXY'PHARM

829 rue Marcel Paul
94500 Champigny-sur-Marne
commercial@oxypharm.net
T : +33 1 45 18 78 70

www.oxypharm.net

